
DIGITAL MICROSCOPE

321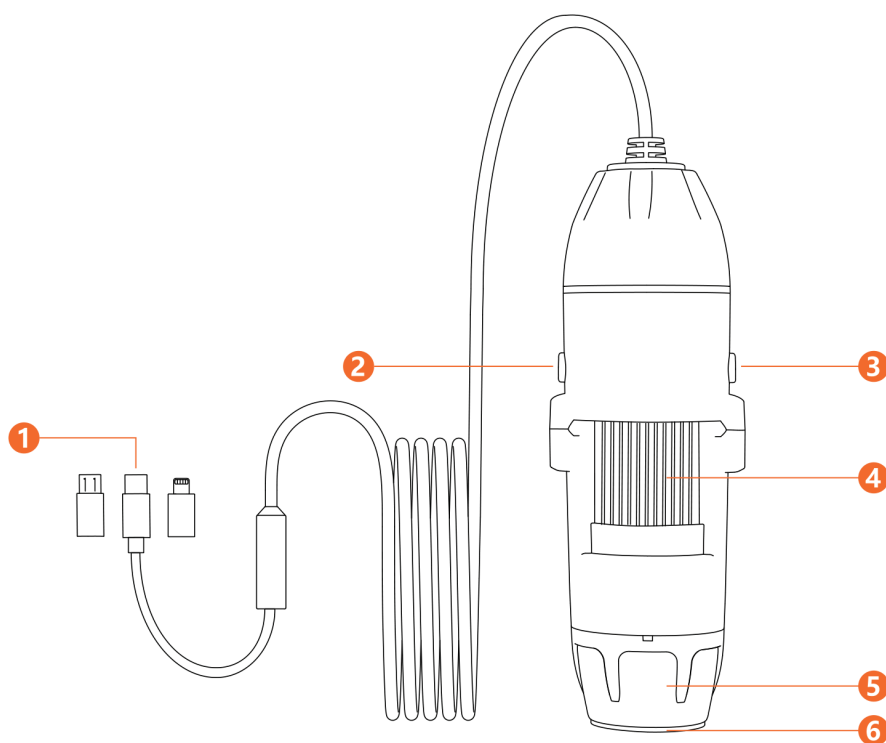

Digital Microscope

1600× magnification · 0–90 mm focus

User Manual

Multilanguage Edition

Čeština · Slovenčina · Magyar
Română · Deutsch · Български
Polski · Slovenščina · English

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321

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
DIGITAL MICROSCOPE

321

Microscope for direct connection to a phone

1600× magnification · 0–90 mm focus · records to your phone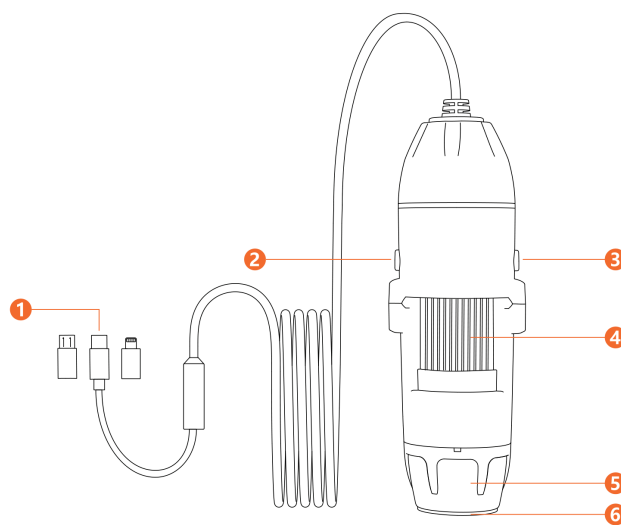

User Manual

Please read this manual before using the device and keep it for future reference.

Contents

1	Important Notice	3
2	Product Introduction	3
3	Product Description	4
4	Installing the app	4
5	Connecting to a phone	5
6	Using the app	6
7	Frequently asked questions	7
8	Specifications	8

1. Important Notice

1. Before first use, remove the protective cover in front of the lens.
2. The device combines a digital and an optical microscope. The actual magnification depends on the captured image.
3. The optimal focal length of the microscope is 0 to 90 mm, and different object distances correspond to different magnifications. To view an object clearly, adjust the focus by rotating the focus wheel.
4. Do not disassemble the device if it is damaged. Contact the place of purchase or your dealer for repair service.
5. For best results, clean the connector on your phone or tablet (Lightning, Type-C or Micro-USB) before use to prevent connection issues.
6. The device is intended to be used under adult supervision.
7. The device is intended for use with smartphones and tablets only. The manufacturer is not liable for damage caused by improper use.


2. Product Introduction

The device is a high-definition microscope that connects directly to the phone connector and displays the image in an app. It takes photographs and records video and works with both Apple phones and phones running Android.

Where the device is used

- textile inspection in the textile industry
- print inspection
- inspection of printed circuit boards and precision mechanisms
- education
- hair inspection
- skin inspection
- observation of micro-organisms
- inspection of jewellery and collectors' coins
- as a visual aid for the partially sighted

3. Product Description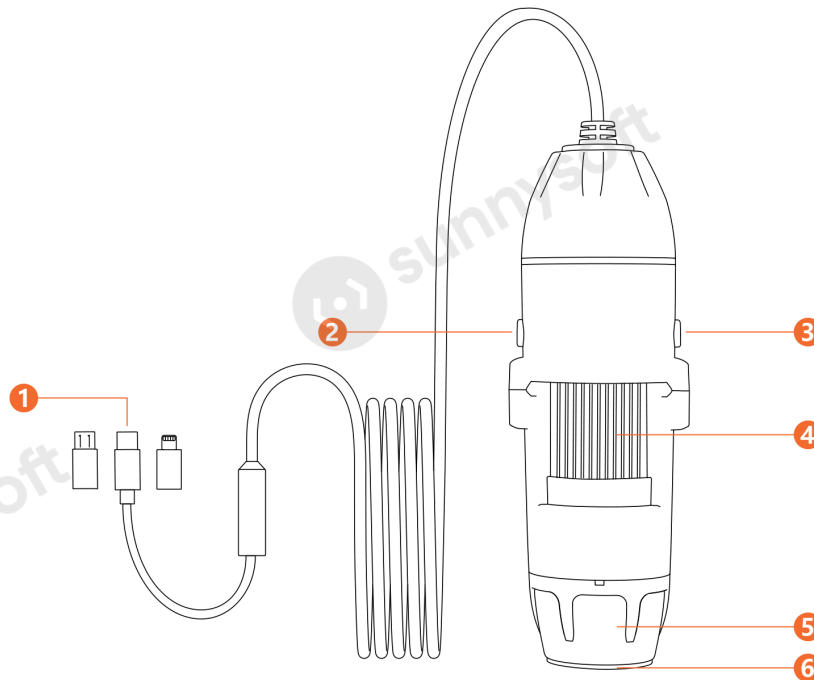

Parts of the device.

- | | |
|----------------------------------|--|
| 1 Micro USB / Type-C / Lightning | 4 Focus wheel |
| 2 Shutter button | 5 Condenser lens |
| 3 Lighting adjustment button | 6 Protective cover (remove before use) |

4. Installing the app

The device works with devices running iOS and with devices running Android that support the OTG function. The app is called "SUP-ANESOK".

Apple phones and tablets (iOS)


Search the App Store for the "SUP-ANESOK" app, then download and install it. iOS 10.0 or later is required.

Phones and tablets running Android

Search Google Play for the "SUP-ANESOK" app, then download and install it. Android 6.0 or later is required.

5. Connecting to a phone

Choose the adapter that matches the connector on your phone and connect the device. The LED on the device lights up. If it does not, check that the connector is fully inserted. On devices running Android, also check in the settings that the OTG function is enabled.

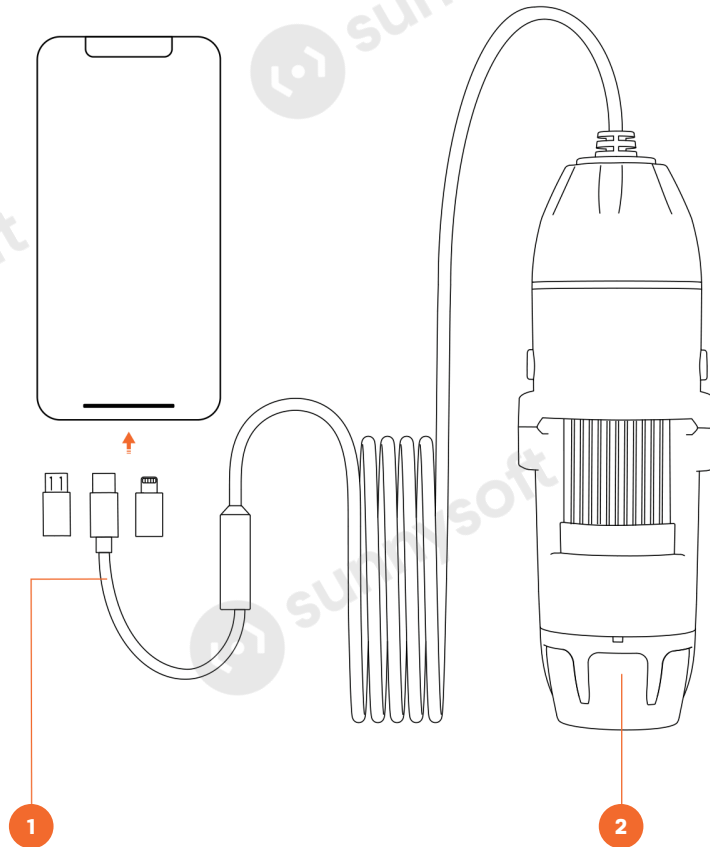
Connecting the device to a phone.

1 Micro USB / Type-C / Lightning

2 LED light location

Removing the protective cover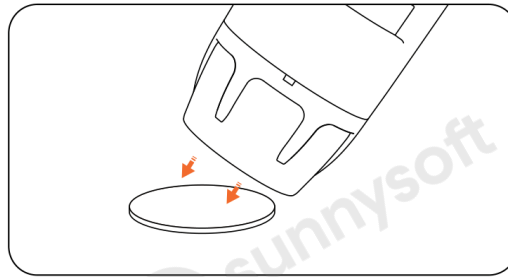

Removing the protective cover from the lens.


Notice

Remove the protective cover before use, otherwise it will obstruct the view.

6. Using the app

Home screen and imaging screen. The app is in English.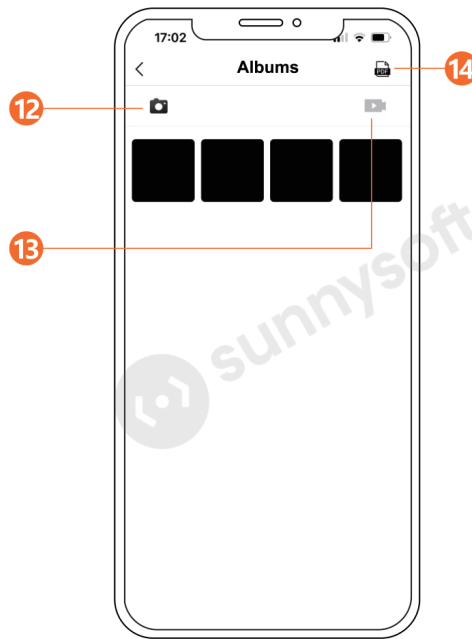

Album of captured photos and videos.

- | | |
|----------------------------------|------------------------------------|
| 1 Settings | 8 Open the album |
| 2 Connection status | 9 Image comparison |
| 3 Go to the imaging screen | 10 Colour and black-and-white mode |
| 4 Open the album | 11 Magnification |
| 5 Back | 12 Browse photos |
| 6 Rotate the image | 13 Browse videos |
| 7 Take a photo or record a video | 14 Create a PDF document |

7. Frequently asked questions

Which systems is the device compatible with?

Android 6.0 and later (with the OTG function) and iOS 10.0 and later.

Why does the app not display an image?

Check whether the LED is on. If it is not, reconnect the device. Also check that the OTG function is enabled in the phone settings. Finally, clean the connector on your phone – see the manual of your phone for instructions.

Why is the image blurred?

Rotate the focus wheel left and right to get the sharpest image and choose a suitable distance from the object. Also make sure the protective cover in front of the lens has been removed.

What causes the static black spots in the image?

Dust, in most cases. Try tapping the device lightly to loosen the dust. If the spots remain, contact your dealer.

Why is the image displayed with a delay?

Make sure the phone battery is charged above 20 %.

8. Specifications

Sensor	2.0 MP
Magnification	up to 1600×
Photo resolution	1920 × 1440
Video resolution	1920 × 1440
Focus	manual, 0 to 90 mm
Photo format	JPG
Video format	MP4
Compatibility	Android 6.0 and later, iOS 10.0 and later

Notice

The photo and video resolution is higher than the resolution of the sensor itself, so the image is interpolated. The values given are taken from the manufacturer's documentation.

Markings on the product

The product carries the CE marking, complies with the RoHS directive on the restriction of hazardous substances and is subject to separate collection of electrical waste – do not dispose of it with unsorted waste. Made in China.

The English version of this manual is a translation of the manufacturer's original manual. The photographs used in the manual are illustrative only and may not correspond exactly to the product.

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